

Pallet Balance Instructions

Use and maintenance methods

1. When using a balance, it should be placed on a flat surface.
2. Move the cursor to the “0” point on the left end of the scale: the pointer should be aligned with the center line of the scale head to achieve balance, otherwise the balance nut on the lever should be rotated to adjust the balance.
3. Place the weights on the right pan and the object to be weighed on the left pan. Try to place the object and the weights in the center of the pan.
4. The balance and weights should be properly stored and should not be placed in places susceptible to moisture and oxidation, and should be kept away from acids, alkalis, greases and oils.
5. The balance and weights should always be kept dry and clean. They must be cleaned with a soft brush before use. The balance and weights should be checked once after use and put back into the box immediately.
6. If the balance or weights are damaged, oxidized, or if there is any doubt about the weight, they should be sent to the manufacturer or repair department for inspection and only used after passing the inspection.

Performance Indicators

Model	Maximum weight (g)	Minimum graduation value (g)	Value scale maximum value (g)
11003	100	0.1	0.1
<i>11004</i>	<i>200</i>	<i>0.2</i>	<i>0.2</i>
11005	500	0.5	0.5
11006	500	1	1
JPT-1	1000	1	1

Note: The bolded and italicized model is the one that came with this kit.